


# TUBE SEALER CR6

**Tube Sealer CR6** is a battery operated sealing unit for EVA, PVC/EVA coextrusion, and PVC blood bag tubes or sets for cryopreservation.

**OriGen<sup>®</sup>**  
BIOMEDICAL

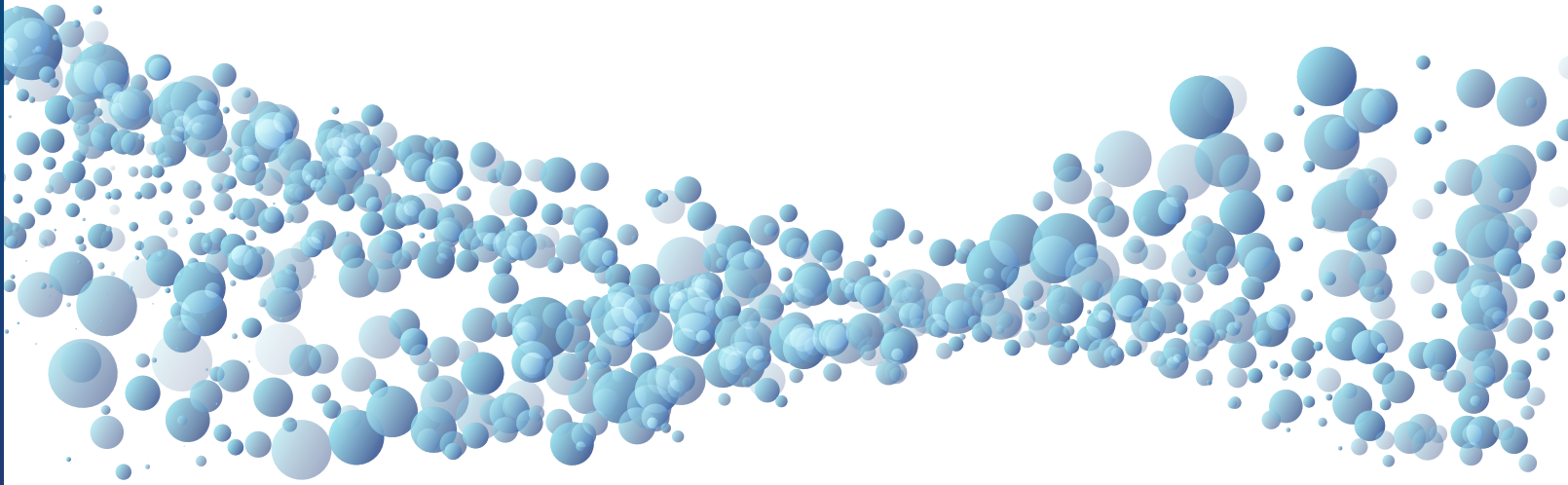

## HIGH QUALITY SEALS OF ORIGEN TUBING

The Tube Sealer CR6 is capable of sealing tubing including OriGen Biomedical<sup>®</sup> EVA, EVA/PVC coextrusions, and PVC tubing with an outer diameter of 6mm. The Tube Sealer CR6 creates reliable, thick seals that can be easily inspected and separated. It can be used with CryoStore Freezing Bags. The device is able to deliver approximately 1,500 seals per charge in PVC tubes.

## PORTABLE

The Tube Sealer CR6 has a power unit with batteries and control unit monitored by microprocessor. The unit comes with a battery charger as well.

## EASY TO USE

Simply place the tube between the sealing handle's electrodes and close the lever. The light on the handle will turn on when sealing begins and will turn off when it is complete. Remove the tube after one additional second. Follow bag recommendations for proper seal placement.

## 1-2 SECOND SEAL TIME

Seal time is an automatic 1-2 second seal time. It is recommended that you conduct less than one seal every three seconds or a maximum of 50 seals in a sequence with a 15 minute rest after that.


Tube Sealer CR6 is comprised of three components:

- Sealing handle with 6' (1.9m) coax cable
- Power unit with batteries and control unit monitored by microprocessor
- Battery charger



# TUBE SEALER CR6

**Tube Sealer CR6** is a battery operated sealing unit for EVA, PVC/EVA coextrusion, and PVC blood bag tubes or sets for cryopreservation.

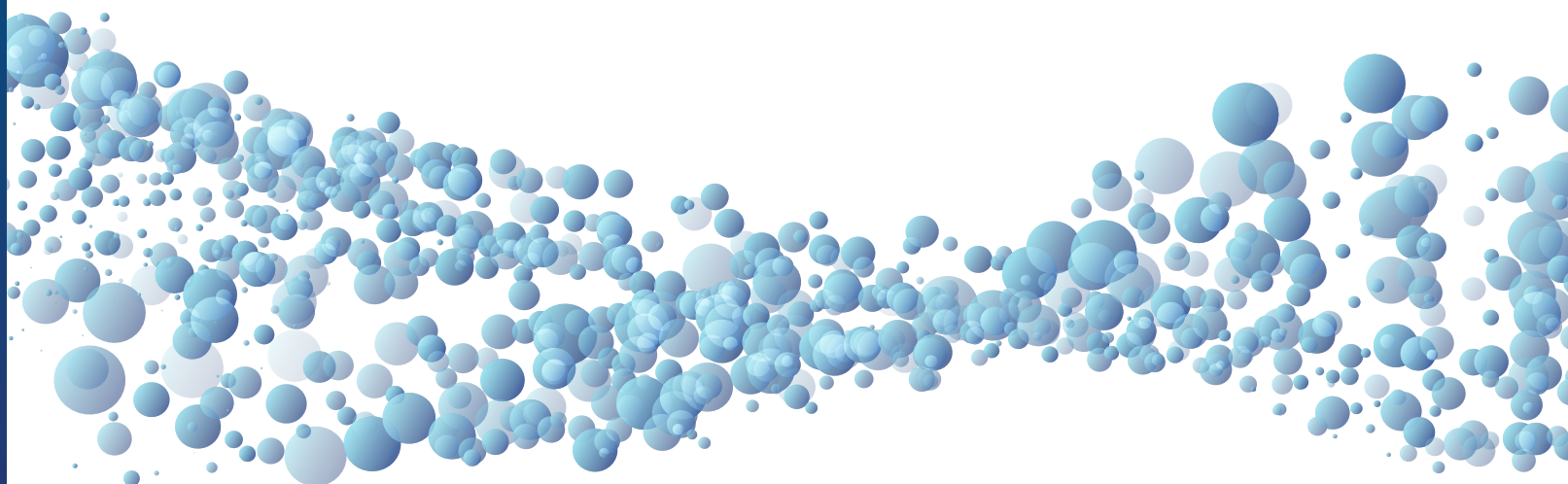

## TECHNICAL DATA

**Batteries:** NiMH battery pack 24V / 3,8Ah


**Seal Time:** Automatic 1-2 Sec

**Dimensions:** 266 x 119 x 88 mm (L x W x H) including HF contact.

**Weight:** 4.7lb (2.14 Kg) including battery

**Operation:** Recommended: Less than 1 seal every three seconds during continuous use or maximum 50 seals in sequence and a 15 minute rest after that.

PART CODE	PRODUCT DESCRIPTION
CR6	CR6 Sealer with smooth seal
CR6P	CR6 Sealer with perforated seam through the seal

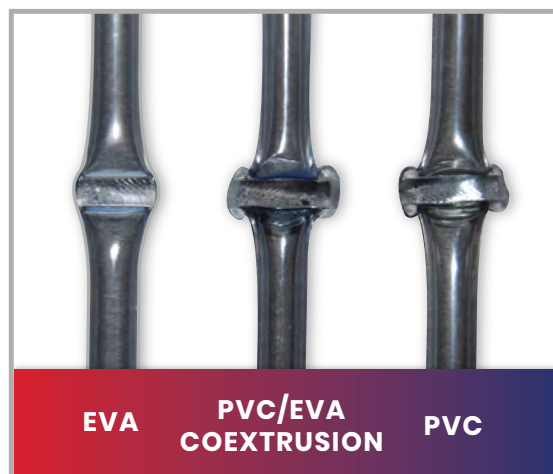

## ORDERING INFORMATION

To place an online order, please fill out [this form](#).

## WARRANTY INFORMATION

Limited Equipment Warranty. Ljungberg & Kögel AB (L&K) hereby guarantees the original buyer that Tube Sealer CR6 is manufactured in a professional and quality manner, and will be free from all faults during a period of one year from the date of delivery from OriGen. L&K will without cost for the customer, repair or replace the equipment that is faulty. The warranty is not valid for the batteries since they are considered to be expendable supplies.

7000 Burluson Rd, Bldg D | Austin, Texas 78744 | 512.474.7278 | Toll Free 800.233.9014 | sales.us@origen.com | origen.com

FDA Registered | ISO 13485 Certified | Manufactured to GMP Guidelines

ML01R05:2023-5

© 2023 OriGen Biomedical. All rights reserved. OriGen Biomedical® is a registered trademark of OriGen Biomedical, Inc